


FieldMaxII-P™

User Manual



**User Manual
FieldMaxII-P
Laser Energy Meter**

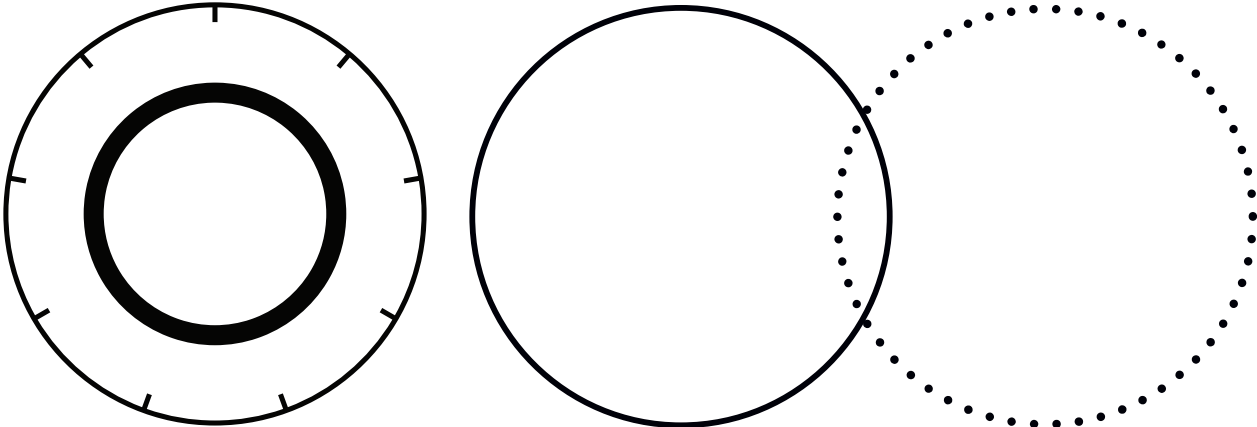

27650 SW 95th Ave.
Wilsonville, OR 97070

This document is copyrighted with all rights reserved. Under the copyright laws, this document may not be copied in whole or in part or reproduced in any other media without the express written permission of Coherent, Inc. Permitted copies must carry the same proprietary and copyright notices as were affixed to the original. This exception does not allow copies to be made for others, whether or not sold, but all the material purchased may be sold, given, or loaned to another person. Under the law, copying includes translation into another language.

Coherent, the Coherent Logo, and FieldMaxII are trademarks or registered trademarks of Coherent, Inc. All other trademarks or registered trademarks are the property of their respective owners.

Patents referenced in this user manual were active as of the printing date of the manual (see last page for date). The patents referenced herein may have expired. You are advised to check to see if the patents are still active: <http://portal.uspto.gov/external/portal/pair>.

Every effort has been made to ensure that the data given in this document is accurate. The information, figures, tables, specifications and schematics contained herein are subject to change without notice. Coherent makes no warranty or representation, either expressed or implied with respect to this document. In no event will Coherent be liable for any direct, indirect, special, incidental or consequential damages resulting from any defects in its documentation.

Technical Support

In the US:

Should you experience any difficulties with your product or need any technical information, please visit our website: www.Coherent.com. Additional support can be obtained by contacting our Technical Support Hotline at 1.800.343.4912 (1.503.454.5700 outside the U.S.), or e-mail LSMupport@Coherent.com.

If you call outside our office hours, your call will be taken by our answering system and will be returned when the office reopens.

If there are technical difficulties with your product that cannot be resolved by support mechanisms outlined above, e-mail, or telephone Coherent Technical Support with a description of the problem and the corrective steps attempted. When communicating with our Technical Support Department via the web or telephone, the Support Engineer responding to your request will require the model and serial number of your system.

Outside the US:

If you are located outside the U.S., visit our website for technical assistance or contact our local service representative. Representative phone numbers and addresses can be found on the Coherent website: www.Coherent.com.

Coherent provides telephone and web technical assistance as a service to its customers and assumes no liability thereby for any injury or damage that may occur contemporaneous with such services. These support services do not affect, under any circumstances, the terms of any warranty agreement between Coherent and the buyer. Operation of any Coherent laser with any of its interlocks defeated is always at the operator's own risk

Table of Contents

1 Introduction	i
1.1 Signal Words and Symbols in this Manual	i
1.1.1 Signal Words	i
1.1.2 Symbols	ii
1.2 Preface	iii
1.3 Safety and Compliance	iii
1.4 Publication Updates	iii
2 Safety and Compliance	1
2.1 Compliance	2
2.1.1 CE Marking	2
2.1.2 Electromagnetic Compatibility	2
2.1.3 Environmental Compliance	3
2.1.4 RoHS Compliance	3
2.1.4.1 China RoHS Compliance	3
2.1.5 Waste Electrical and Electronic Equipment (WEEE, 2002)	4
2.1.6 Battery Directive	4
2.2 Export Control Laws Compliance	4
2.3 Declaration of Conformity	5
3 Description	7
3.1 Features	7
3.2 Front Panel	8
3.2.1 Buttons	9
3.2.2 Display	10
3.3 Right Side Panel	11
3.3.1 USB Connector	11
3.3.2 Sensor Connector	12
3.4 Left Side Panel	12
3.4.1 Analog Out Connector	12
3.4.2 Power Jack Connector	12
3.5 AC Adapter	12
4 Quick Setup and Getting Started	15
4.1	
Install Software and Drivers	15
4.1.1 Install Software	15
4.1.2 USB Driver Installation	16
4.2 Measure Power With a Thermopile or Optical Sensor	16
5 Operation	17
5.1 Tuning Mode	17
5.1.1 Tuning Needles	17
5.1.2 Zone Indicator Bars	18

5.2 Annunciators	20
5.2.1 Temperature	20
5.2.2 AUTO	20
5.2.3 Range Hint	20
5.2.4 AVG	20
5.2.5 ATTEN	20
5.2.6 RMT	21
5.2.7 Battery	21
5.2.8 Hold	21
5.3 Fault Display Indicators	21
5.4 Invalid and not Available Data Conditions	22
5.4.1 Invalid Data	22
5.4.2 Not Available Data	23
5.5 Setup Parameters	23
5.5.1 Setup/Local	23
5.5.2 Stat Button	24
5.5.3 Wave Button	24
5.5.4 Area Button	25
5.5.4.1 Area Correction and Zeroing	25
5.5.5 Avg Button	26
5.5.6 Atten Button	26
5.5.7 Hold	27
5.6 Button Functions	27
5.6.1 Auto	27
5.6.2 Zero Button	27
5.6.3 Power Switch and Backlight Toggle Button	28
5.6.4 Up and Down Arrows	28
5.6.5 Left and Right Arrows	28
5.7 Statistics Mode	28
5.7.1 Invalid Data	29
5.8 Measurement Display and Range Selection	29
5.8.1 Manual Ranging Mode	31
5.8.2 Auto vs. Manual Ranging Mode	31
5.8.3 Zero the Sensor	31
5.8.4 Negative Power Display	32
6 Host Interface	33
6.1 Overview	33
7 Calibration and Warranty	35
7.1 Calibration	35
7.1.1 Scope of Calibration	35
7.1.2 Re-certify Once Per Year	35
7.1.3 Coherent Calibration Facilities and Capabilities: ISO 17025 Accredited	36
7.2 Limited Warranty	36
7.3 Warranty Limitations	36
8 Extended Warranty Program	39

Appendix A: Maintenance	41
A.1 Battery Replacement.....	41
Appendix A: Service & Support	43
A.1 Technical Support.....	43
A.1.1 Support in the USA and North America.....	43
A.1.2 International Support.....	44
A.2 Obtain Service.....	44
A.3 Product Shipment Instructions	45
Appendix B: Specifications	47
Index	49

List of Figures

2-1.	China RoHS Label.....	3
2-2.	WEEE Label.....	4
3-1.	FieldMax II TO.....	7
3-2.	Front Panel.....	8
3-3.	LCD Display.....	10
3-4.	Right Side Panel.....	11
3-5.	Left Side Panel.....	12
4-1.	Measure Power With a Thermopile or Optical Sensor.....	16
5-1.	Location of Tuning Needles and Zone Indicator Bars.....	18
5-2.	Current Scale Mid-Range.....	18
5-3.	Zone Indicator Bars.....	18
5-4.	Comparison of Zone Indicator Bars to Full Scale Measurement.....	19
5-5.	Tuning Mode Example - Full Scale.....	19
5-6.	Auto Ranging.....	31
A-1.	Optional Alkaline Battery Replacement.....	42

List of Tables

5-1.	Fault Codes.....	23
5-2.	Full Scale Range Settings.....	30
A-1.	Batteries Contained in this Product.....	43
B-1.	Specifications.....	49

1 Introduction

1.1 Signal Words and Symbols in this Manual

This documentation may contain sections in which particular hazards are defined or special attention is drawn to particular conditions. These sections are indicated with signal words in accordance with ANSI Z-535.6 and safety symbols (pictorial hazard alerts) in accordance with ANSI Z-535.3 and ISO 7010.

1.1.1 Signal Words

Four signal words are used in this documentation: **DANGER**, **WARNING**, **CAUTION** and **NOTICE**.

The signal words **DANGER**, **WARNING** and **CAUTION** designate the degree or level of hazard when there is the risk of injury:

DANGER!

Indicates a hazardous situation that, if not avoided, will result in death or serious injury. This signal word is to be limited to the most extreme situations.

WARNING!

Indicates a hazardous situation that, if not avoided, could result in death or serious injury.

CAUTION!

Indicates a hazardous situation that, if not avoided, could result in minor or moderate injury.

The signal word '**NOTICE**' is used when there is the risk of property damage:

NOTICE

Indicates information considered important, but not hazard-related.

Messages relating to hazards that could result in both personal injury and property damage are considered safety messages and not property damage messages.

1.1.2

Symbols

The signal words **DANGER**, **WARNING**, and **CAUTION** are always emphasized with a safety symbol that indicates a special hazard, regardless of the hazard level:



This symbol is intended to alert the operator to the presence of important operating and maintenance instructions.



This symbol is intended to alert the operator to the danger of exposure to hazardous visible and invisible laser radiation.



This symbol is intended to alert the operator to the presence of dangerous voltages within the product enclosure that may be of sufficient magnitude to constitute a risk of electric shock.



This symbol is intended to alert the operator to the danger of Electro-Static Discharge (ESD) susceptibility.



This symbol is intended to alert the operator to the danger of crushing injury.



This symbol is intended to alert the operator to the danger of a lifting hazard.

1.2

Preface

This manual contains user information for the FieldMaxII-P™ laser energy meter.

1.3

Safety and Compliance

Read and obey all safety and compliance information provided in the next chapter before equipment is set up and operated.

1.4

Publication Updates

To view information that may have been added or changed since this publication went to print, connect to www.Coherent.com/resources.

2 Safety and Compliance

Carefully review the following safety information to avoid personal injury and to prevent damage to this meter or any sensor connected to it. Except for replaceable batteries—discussed under 'AC Adapter' (p. 12)—there are no user-serviceable parts in the FieldMaxII-P meter. For service information, refer to 'THIS WARRANTY IS EXCLUSIVE IN LIEU OF ALL OTHER WARRANTIES WHETHER WRITTEN, ORAL, OR IMPLIED. COHERENT SPECIFICALLY DISCLAIMS THE IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE. IN NO EVENT SHALL THE COMPANY BE LIABLE FOR ANY INDIRECT, INCIDENTAL OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH ITS PRODUCTS.' (p. 39).



WARNING!

Use only the power cord specified for the meter. The grounding conductor of the cord must be connected to earth ground.



WARNING!

Do not operate the meter if its panels are removed or any of the interior circuitry is exposed.



WARNING!

Do not operate the meter in wet or damp conditions, or in an explosive atmosphere.



NOTICE

Operate the meter only within the specified voltage range.



NOTICE

Do not apply a voltage outside the specified range of the input connections.



NOTICE

Do not operate the meter if there are suspected failures. Refer damaged units to qualified Coherent service personnel.

2.1 Compliance

This section describes compliance with various government requirements for safety, environmental regulations, and control law.

For additional information about Coherent Compliance, please visit:

<https://www.coherent.com/company/environmental>

2.1.1 CE Marking

The European Community requirements for product safety are specified in the Low-Voltage Directive (LVD) (published in 2014/35/EU).

This Directive requires that lasers comply with the standard EN 61010-1/IEC 61010-1 “Safety Requirements For Electrical Equipment For Measurement, Control and Laboratory Use” and EN 60825-1/IEC 60825-1 “Safety of Laser Products”. Compliance with the European requirements is certified by CE Marking.

2.1.2 Electromagnetic Compatibility

The primary issue for electromagnetic compatibility is to design covers, shielding, grounding, routing of electrical cable assemblies, and control elements with the proper safety features for a complete system.

The European requirements for Electromagnetic Compliance (EMC) are specified in the EMC Directive (published in 2014/30/EU).

Class A:

Conformance (EMC) concerning emission and immunity is achieved through compliance with the harmonized standard EN 61326-1_Ed2:2013 (IEC 61326-1_Ed2:2012 for Electrical Requirement for Measurement, Control and Laboratory for Class A.

Class B:

Conformance to the EMC requirements is achieved through compliance with the harmonized standards EN55011:2009 for emission and EN61000-6-2:2006 for immunity.

2.1.3 Environmental Compliance

This section describes compliance with various environmental regulatory directives to identify hazardous substances.

2.1.4 RoHS Compliance


The European Union RoHS Directive EN 50581:2012 restricts the use of certain hazardous substances in electrical and electronic equipment.

Coherent product(s) conform to all applicable requirements of the EU-RoHS Directive (2011/65/EU) and Amendment Directive (EU) 2015/863. Compliance Declarations are available upon request.

2.1.4.1 China RoHS Compliance

The China RoHS (Restriction of Hazardous Substances) Regulation SJ/T 11364-2014 restricts the use of certain hazardous substances in electrical and electronic equipment. The China RoHS Regulation applies to the production, sale, and import of products into the Peoples Republic of China.

Any hazardous substances in Coherent products (if applicable) are listed on the product label, as shown in the example shown in Figure 2-1:

产品中有害物质的名称及含量							
部件名称 Part Name	有害物质 Hazardous Substances						
	铅 (Pb)	汞 (Hg)	镉 (Cd)	六价铬 (Cr(VI))	多溴联苯 (PBB)	多溴二苯醚 (PBDE)	
印刷电路板组装 Printed Circuit Board Assembly	X	O	O	O	O	O	
电源 Power Supply	X	O	O	O	O	O	
电源线 Power Cord	X	O	O	O	O	O	

本表格依据 SJ/T 11364 的规定编制
 O: 表示该有害物质在该部件所有均质材料中的含量均在 GB/T 26572 规定的限量要求以下。
 X: 表示该有害物质至少在该部件的某一均质材料中的含量超出 GB/T 26572 规定的限量要求。

Figure 2-1. China RoHS Label

The China RoHS Regulation also requires that the date of manufacture be identified. This information is provided on the *Certificate of Calibration* shipped with each product.

2.1.5 Waste Electrical and Electronic Equipment (WEEE, 2002)

Coherent product(s) conform to all applicable requirements of the EU Waste Electrical and Electronic Equipment (WEEE)- Directive (2012/19/EU). WEEE management also covers EU Directive 2006/66/EC-EU Battery Directive and Directive 94/62/EC on Packaging and Packaging Waste. Do not dispose of these products or packaging as unsorted municipal waste.

Coherent joins approved compliance organizations to meet its collection and recycling obligations. For further information, please contact:

Email: info@rene-europe.com
Phone: +49 (0) 8266-869806
Website: www.rene-europe.com

This directive is represented by a crossed-out garbage container label, shown in Figure 2-2.

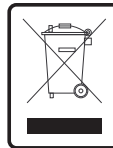


Figure 2-2. WEEE Label

2.1.6 Battery Directive

The batteries used in this product are in compliance with the EU Directive 2006/66/EC (“EU Battery Directive”). For a list of batteries used with the product, refer to Table A-1, ‘Batteries Contained in this Product,’ (p. 43).

2.2 Export Control Laws Compliance

It is the policy of Coherent to comply strictly with U.S. export control laws.

Export and re-export of lasers manufactured by Coherent are subject to U.S. Export Administration Regulations, which are administered by the Commerce Department. In addition, shipments of certain components are regulated by the State Department under the International Traffic in Arms Regulations.

The applicable restrictions vary depending on the specific product involved and its destination. In some cases, U.S. law requires that U.S. Government approval be obtained prior to resale, export or re-export of certain articles. When there is uncertainty about the obligations imposed by U.S. law, clarification must be obtained from Coherent or an appropriate U.S. Government agency.

Products manufactured in the European Union, Singapore, Malaysia, Thailand: These commodities, technology, or software are subject to local export regulations and local laws. Diversion contrary to local law is prohibited. The use, sale, re-export, or re-transfer directly or indirectly in any prohibited activities are strictly prohibited.

2.3 Declaration of Conformity

Declaration of Conformity certificates are available upon request.

3 Description

Thank you for purchasing the FieldMaxII-P™—a versatile, easy-to-use digital energy meter designed for field service and production applications.



Figure 3-1. FieldMax II P

3.1 Features

Below is a list of specific features included in the FieldMaxII-P meter:

- 73 x 58 mm backlit LCD display
- Fast and effective laser tuning mode
- Works with pyroelectric sensors
- Measures energy up to 300 pps
- Intuitive soft key-driven user interface
- USB 1.1 interface
- Portable AC/DC operation
- Compact, rugged enclosure with stand

The versatile FieldMaxII-P measures:

- Energy: J, J/cm²
- Frequency: Hz
- Full statistics: max, min, mean, and standard deviation

3.2 Front Panel

The front panel—Figure 3-1, below—includes a liquid crystal display (LCD) and buttons that are used to enter parameters, select modes, and change ranges.

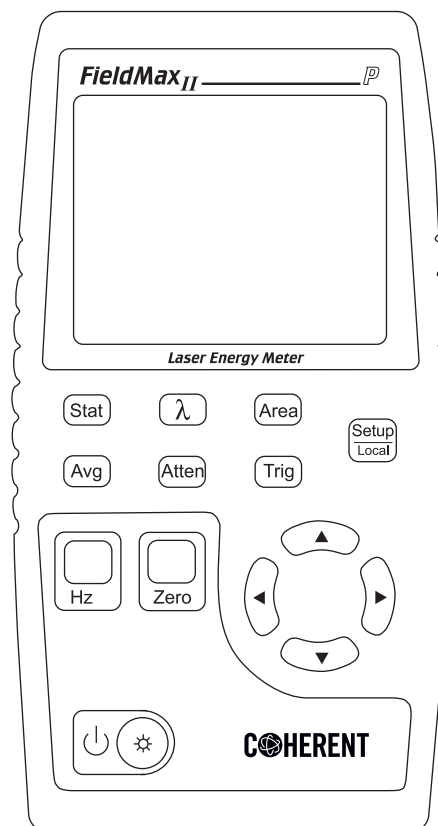



Figure 3-2. Front Panel

3.2.1

Buttons

The following buttons are on the front panel of the FieldMaxII-P:

- Setup/Local—starts or ends a parameter edit cycle. This button is also used to cancel the front panel lockout when Remote Control is active.
- Stat—statistics processing parameter
- Wave (λ)—wavelength compensation parameter
- Area—area correction parameter
- Avg—display smoothing parameter
- Atten—attenuation correction parameter
- Trig—trigger level parameter
- Hz—Rep Rate Display mode
- Zero—start batch. When Statistics mode is not active, this button can also be used to restart a batch.
- —Power Switch/Backlight Toggle button
- Up Arrow (\blacktriangle)—field adjust or range select
- Down Arrow (\blacktriangledown)—field adjust or range select
- Left Arrow (\blacktriangleleft)—field select
- Right Arrow (\blacktriangleright)—field select

For detailed information about each of these buttons, refer to 'Operation' (p. 19).

3.2.2 Display

The LCD display provides visual measurement information. Figure 3-3 shows all the possible segments that may appear on the display.

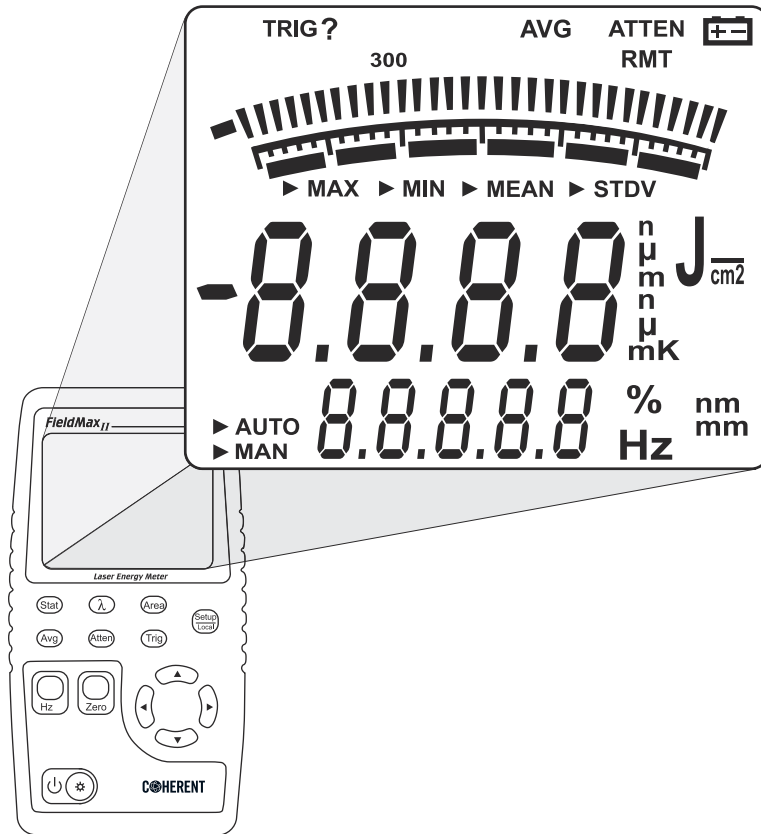


Figure 3-3. LCD Display

NOTICE

The type of sensor being used and the individually-selected settings determine what type of information will actually appear on the display

Information that appears on the display is divided into the groups described in the following list. Figure 3-1 shows the general location of each group).

- Annunciators: TRIG, Range Hint, AVG, ATTEN, RMT, and Battery
- Digital tuning feature

- Tuning meter scale
- Statistical parameters: MAX, MIN, MEAN, and STDV
- Numeric measurement value—large numeric characters
- Measurement units and engineering prefixes
- Numeric data entry, batch count, parameter settings, and Hertz measurement values
- Data entry units, Hertz units, and current parameter units

For detailed information about these settings, refer to 'Button Functions' (p. 28).

3.3 Right Side Panel

The right side panel contains the USB and Sensor connectors—refer to Figure 3-4.

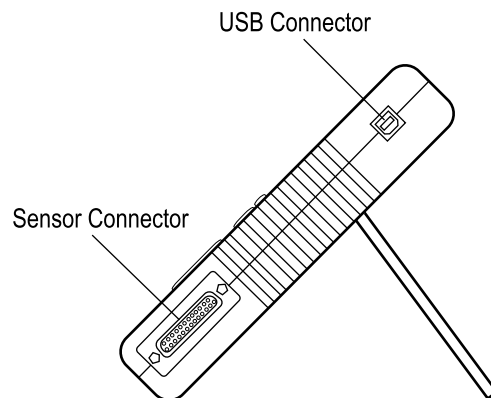


Figure 3-4. Right Side Panel

3.3.1 USB Connector

Attachment of the supplied cable to this standard USB connector allows communication between FieldMaxII-P and a computer with a USB interface.

3.3.2 Sensor Connector

Use this connection to attach a DB-25 SmartProbe connector or adapter.

3.4 Left Side Panel

The left side panel contains the Analog Out and Power Jack connectors. Refer to Figure 3-5.

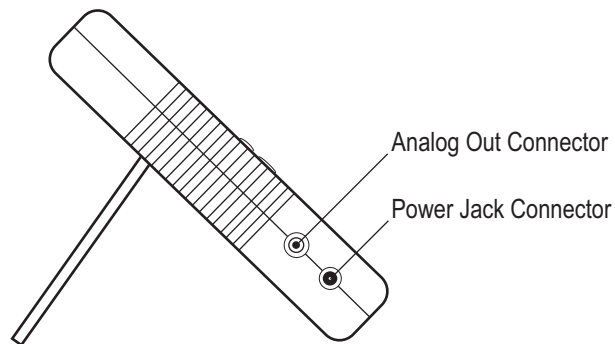


Figure 3-5. Left Side Panel

3.4.1 Analog Out Connector

When power is set to ON, the Analog Out connector outputs a voltage proportional to the current laser measurement. The output voltage is zero (0) volts when the measured energy is zero (0) or less. The output voltage is the full-scale output voltage when the measured energy is full-scale or over-ranged. The full-scale output voltage (1V, 2V, or 5V) is selected via the host interface. Factory default full-scale output voltage is 2V.

3.4.2 Power Jack Connector

The supplied power cord is connected to this jack.

3.5 AC Adapter

Use of an AC adapter prolongs battery life. FieldMaxII-P automatically senses when an adapter is used.

Description

When batteries rather than an adapter, are used, the battery annunciator flashes if the battery charge is low.

While in battery operation, if a sensor is not connected to the meter, power is automatically set to OFF after ten minutes.

4 Quick Setup and Getting Started

This section presents a 'mini-tutorial' that explains how to connect a sensor to a FieldMaxII-P meter and begin to take measurements within minutes. For in-depth information about product functions for the procedure introduced in this section, refer to 'Operation' (p. 19).

In this section:

- 'Install Software and Drivers' (p. 15)
- 'Measure Energy With a Pyroelectric Sensor' (p. 16)



WARNING!

Follow all laser safety procedures. The laser must be blocked or switched OFF before beginning the procedure described in this section.



WARNING!

Power to the FieldMaxII-P instrument must be OFF before beginning the procedure described in this section.

4.1

Install Software and Drivers

4.1.1

Install Software

For complete software installation instructions, refer to the *FieldMax II Software Installation and Quick Start Guide* (PN 1176436), available on the Coherent website.

If that document is unavailable, go to www.coherent.com/resources and download the Fieldmax II Setup software.

4.1.2 USB Driver Installation

When first connecting the meter to a PC with the USB cable, users are prompted through an installation process. USB drivers are automatically installed onto the computer.

Insert the supplied USB drive when prompted.

NOTICE

Do not exceed the energy density limits of the sensor.

4.2 Measure Energy With a Pyroelectric Sensor

The following figure outlines how to set up a pyroelectric sensor to take an energy measurement. Special Topics

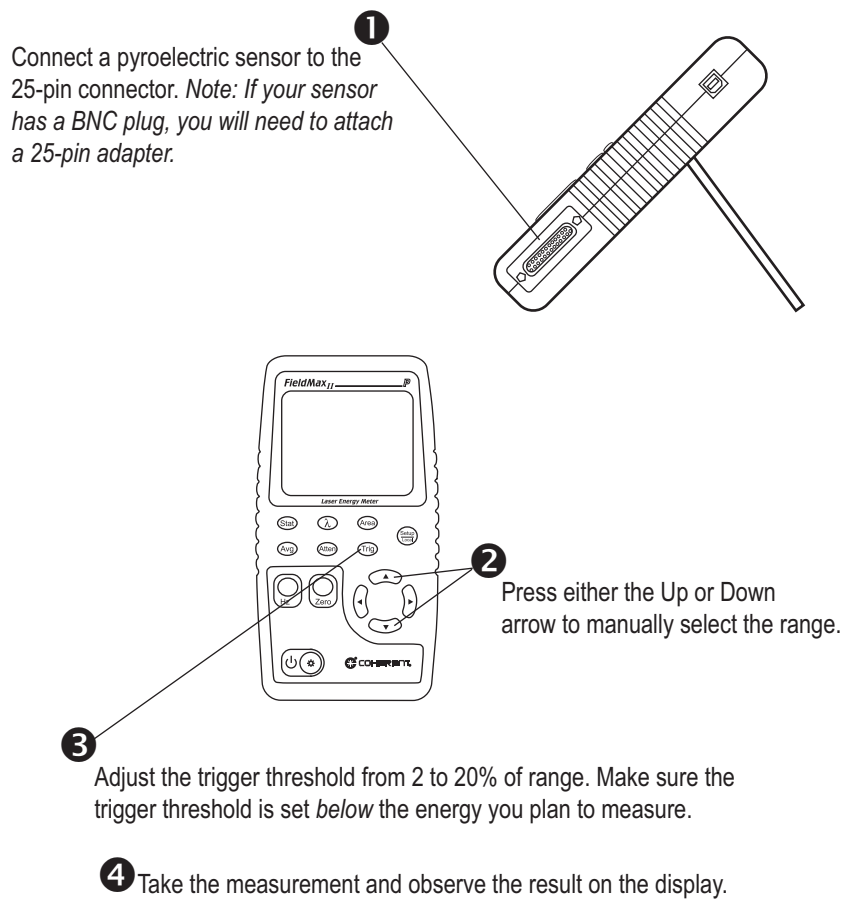


Figure 4-1. Measuring Energy With a Pyroelectric Sensor

This section gives more information on trigger states and modes of operation.

4.2.1 Trigger States and the Trigger Annunciator

Two segments are dedicated to the indication of the trigger state of the meter. These are the TRIG and the '?' segment that follows it.

There are two possible trigger states: 'triggered' and 'wait for trigger.' The triggered state indicates that the meter is capturing laser pulses. The wait for trigger state indicates an absence of pulse triggers. When a pulse is captured, the trigger state becomes active and the wait for trigger state becomes inactive. If no pulses are captured within 1.67 seconds since the last pulse was captured, the wait for trigger state becomes active and the triggered state becomes inactive. The TRIG annunciator reads "TRIG" in the triggered state and "TRIG?" in the wait for trigger state.

5 Operation

This section gives detailed information about control functions and how they are used operate the Fieldmax II:

- 'Tuning Mode' (p. 19)
- 'Annunciators' (p. 22)
- 'Fault Display Indicators' (p. 23)
- 'Invalid and not Available Data Conditions' (p. 24)
- 'Setup Parameters' (p. 24)
- 'Button Functions' (p. 28)
- 'Statistics Mode' (p. 29)
- 'Measurement Display and Range Selection' (p. 29)

5.1 Tuning Mode

As shown in Figure 5-1, below, tuning is visually displayed on the LCD using tuning needles and zone indicator bars.

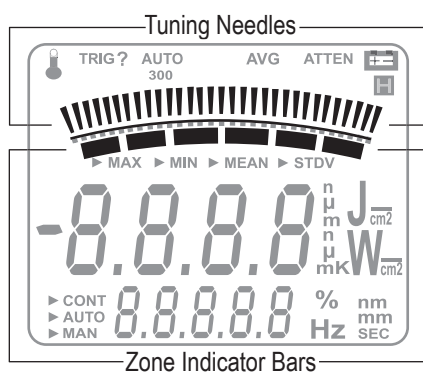


Figure 5-1. Location of Tuning Needles and Zone Indicator Bars

5.1.1 Tuning Needles

Tuning needles—which divide a given tuning zone into thirty ‘increments’—are used to bring laser output to a peak. As the top or bottom of a zone is reached, the tuning needles automatically move to the center of the next zone (see Figure 5-2). Zone indicator bars shows when this happens. Refer to ‘Zone Indicator Bars’ (p. 20) for more information.

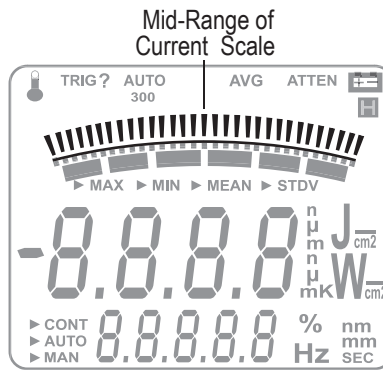


Figure 5-2. Current Scale Mid-Range

5.1.2 Zone Indicator Bars

Zone indicator bars are a series of six segments, as shown in Figure 5-3, below:

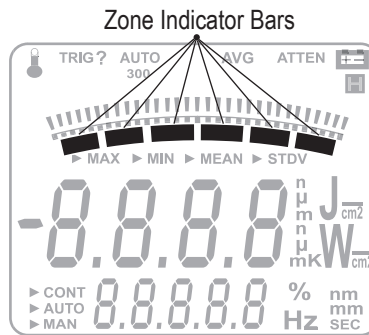


Figure 5-3. Zone Indicator Bars

These bars act as visual indicators while the tuning needles automatically move through zones, and also provide a relative indication of where the measurement falls within the active range.

Zone indicator bars always appear in pairs, with each overlapping zone representing 1/3 of full scale. Figure 5-4 shows how the five zone indicator bars correlate to full scale measurement on the tuning meter scale.

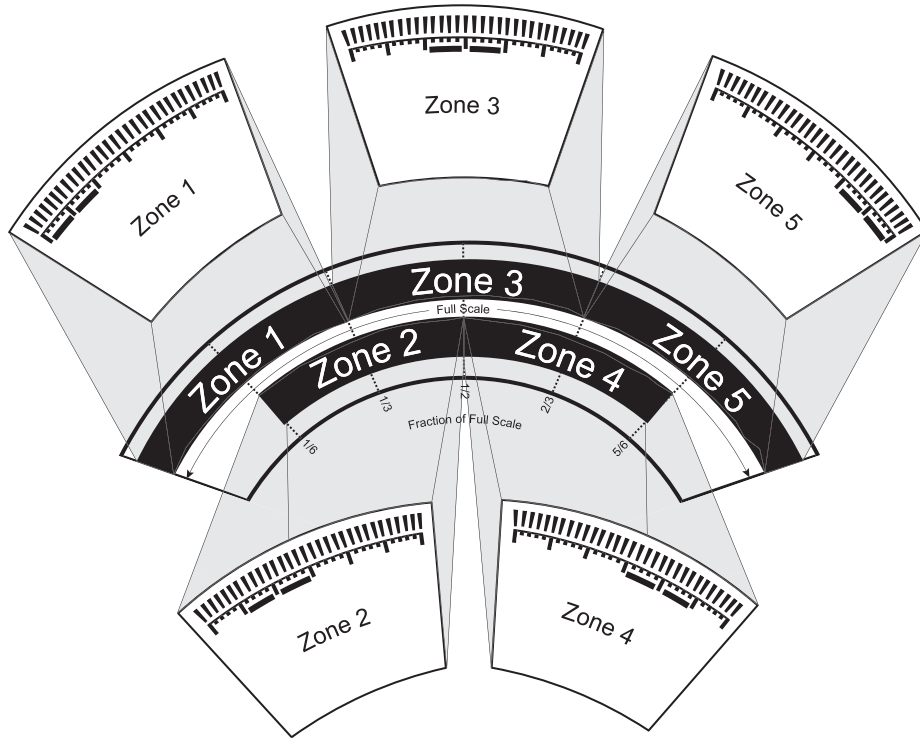


Figure 5-4. Comparison of Zone Indicator Bars to Full Scale Measurement

Figure 5-5 shows an example of how the zone indicator bars overlap on a 30-watt scale:

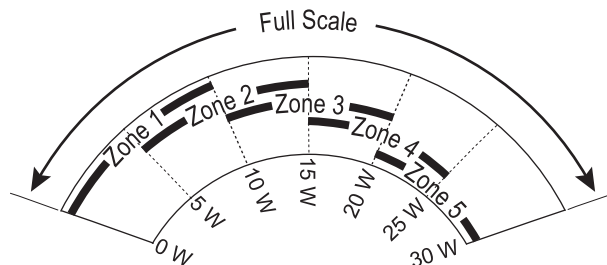


Figure 5-5. Tuning Mode Example - Full Scale

5.2 Annunciators

Annunciators refers to the icon-type symbols that appear on the LCD. Figure 3-3 (p. 10) shows all the annunciators on the FieldMaxII-P meter. The update rate for all annunciators is 3 times per second.

5.2.1 TRIG

Whenever a trigger is detected, “TRIG” displays in the TRIG annunciator position. “TRIG?” displays in the absence of a trigger.

5.2.2 Range Hint

The Range Hint annunciator—towards the top of the LCD, just above the tuning needles—displays ‘3,’ ‘30,’ or ‘300’. These numbers represent the full-scale range currently selected by the user. Range Hint is discussed in more detail under ‘Measurement Display and Range Selection’ (p. 29).

5.2.3 AVG

When AVG (display smoothing) is active, display values are averaged by samples. “AVG” displays when averaging is active. Nothing displays in this position if averaging is not active.


5.2.4 ATTEN

The ATTEN annunciator indicates if attenuation correction is applied to the measurement value. ‘ATTEN’ displays whenever attenuation is active.

5.2.5 RMT

The RMT annunciator indicates that the FieldMaxII meter is currently in Remote Control mode. Remote Control is discussed under ‘Setup/Local’ (p. 24).

5.2.6

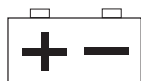Battery

The Battery annunciator flashes whenever the batteries need to be replaced.

The Battery annunciator only appears when the meter is running on battery power, not the AC adapter.

5.3

Fault Display Indicators

FieldMaxII-P is capable of detecting internal and user-induced faults. When a fault is detected, the letters 'Er', followed by a numeric fault code, appear on the display—refer to Table 5-1, below. A fault can be dismissed by pressing any button, or by correcting the cause of the fault.

Table 5-1. Fault Codes

Error condition	Fault code
Unrecognized sensor	1
Sensor communication failure	2
Sensor error	3
Sensor error	4
Sensor/firmware version mismatch (sensor format version exceeds capability of the instrument firmware—firmware upgrade needed)	5
Sample rate fault	6
Hardware fault (detectable hardware error)	20
Data overflow (result of an arithmetic operation that is greater than can be held in the allocated storage)	41
Wrong type of sensor is attached to the instrument	42

As an example, 'Er 4' appears on the display if there is a sensor error. The fault can be dismissed by removing the sensor from the meter, or by pressing any button.

Attaching an unrecognized sensor to the meter creates a special fault condition. This condition is characterized by displaying a sensor fault (1 through 5).

5.4 Invalid and not Available Data Conditions

The update rate for invalid or not available data conditions is three times per second.

5.4.1 Invalid Data

Invalid data is obtained whenever the meter over-ranges. If "OL" appears, start a new batch by pressing the Zero button. All data used to generate a batch result must be valid.

5.4.2 Not Available Data

Measurement data may be unavailable at certain times during meter operation. When data is not available, a series of dashes appears in the measurement area of the display. The following conditions will generate unavailable data:

- Meter is powered on and no pulse triggers are detected.
- Statistics mode is entered and batch data has not been compiled.

5.5 Setup Parameters

This section explains how to select and set user-definable parameters.

5.5.1 Setup/Local

The Setup/Local button serves several purposes:

1. If an edit cycle is not in progress, pressing the Setup/Local button initiates an edit cycle. If the next button pressed represents an edit

parameter (STAT, WAVE, AREA, ATTEN, TRIG, or AVG button), Edit mode is entered and the parameter for the applicable button may be edited using the edit buttons. Additional presses of the Setup/Local button—without first pressing an edit parameter button—are ignored.

2. If a parameter has not been selected, pressing the Setup/Local button cancels an edit cycle.
3. If Edit mode is active and a parameter has been selected, pressing the Setup/Local button commits the adjusted edit parameter to the instrument.
4. Setup/Local is used to cancel the front panel lockout when Remote Control is active.

NOTICE

Remote Control is active when the instrument is connected to a host computer via a USB connection and is communicating with a host application program. While the instrument is in Remote Control mode, all front panel buttons—except Power/Backlight and Setup/Local—are disabled. Pressing the Setup/Local button while in Remote Control mode cancels Remote Control and returns the instrument to Local (all instrument functions available) mode.

The Setup/Local button is also used to edit user-defined parameters. Pressing this button initiates a parameter edit cycle.

A parameter edit cycle consists of:

- Pressing the Setup/Local button to begin the cycle.
- Pressing the parameter button (Stat, Wave (λ), Area, Avg, Atten, or Trig) that needs to be edited. If an edit cycle has been initiated, successive presses of the same edit parameter button will be ignored.
- Using the arrow buttons to select the appropriate field and adjust the data value.
- Pressing the Setup/Local button a second time to end the cycle and commit the new data value.

5.5.2

Stat Button

A press of the Stat button will:

- Enter Statistics mode—if Statistics mode is not active and the button is pressed for less than two seconds.

- Exit Statistics mode.
- Enter Edit mode and select the statistics parameters to be edited (Batch Size and Restart mode) after edit cycle initiation (if Setup/Local is pressed beforehand). Batch size is 2 to 99,999 pulses.
- View the statistics parameters if the button is pressed for two seconds or more.

5.5.3 Wave Button

λ Users can configure FieldMaxII-P to automatically account for any difference between the laser wavelength and the calibration wavelength.

After pressing Setup/Local, the Wave button is used to enter Edit mode and set the wavelength. If an edit cycle has not been initiated, pressing the Wave button will display the wavelength value. The available wavelength range is 1.00 to 99,999 nm. The actual range is sensor-dependent.

NOTICE

If Wavelength compensation information is not programmed into the sensor, users will not be able to change the wavelength data value.

5.5.4 Area Button

This mode allows the measurement of laser energy in terms of fluence.

The parameter is entered as a diameter and assumes a circular beam or aperture. The range for Area mode is 0.01 to 999.99 mm.

Pressing the Area button will:

- Toggle the state of Area Correction mode, if the button is pressed for less than one second. Enter Edit mode and select the beam diameter, if Setup/Local is pressed beforehand.
- View the beam diameter, if the button is pressed for one second or more.

5.5.5 Avg Button

Average mode enables display smoothing, which suppresses variations in the display reading that can make it difficult to read.

NOTICE

Display values are averaged either by time (for thermopile or optical sensors), or by pulses (for pyroelectric sensors).

NOTICE

Display values are averaged by pulses.

The FieldMaxII-P uses a boxcar averaging method, with a boxcar length of 32 pulses.

Pressing the Avg button toggles the state of Average mode.

5.5.6 Atten Button

Attenuation mode allows true measurements with use of an attenuator that has a known attenuation factor. When Attenuation Correction mode is enabled, the measured value is adjusted to indicate the measurement at the attenuator and not the sensor. The range for this mode is 0.01 to 999.99.

Here's an example of how to determine the attenuation correction factor that needs to be set in the FieldMaxII-P meter: If a 1 J laser beam is focused through an attenuator that has an attenuation factor of 50%, then, to get a true laser measurement value, the correction factor in the FieldMaxII-P instrument needs to be set to 2. In other words, since only half the energy of the beam is transmitted through the attenuator, the measured result must be doubled to obtain a true laser measurement.

The Atten button is used to initiate several activities:

- Toggle the state of Attenuation Correction mode, if the button is pressed for less than one second.
- Enter Edit mode and select the attenuation factor, if Setup/Local is pressed beforehand.
- View the attenuation factor, if the button is pressed for one second or more.

5.5.7 Trig Button

After pressing the Setup/Local button, the Trig button is used to enter Edit mode and select the trigger level parameter that will be edited. If an edit cycle has not been initiated, pressing the Trig button will display the trigger level parameter. Trigger has a range of 2 to 20% of full scale. Refer to 'Internal Triggering Mode' (p. 31) for details of the trigger function

5.6 Button Functions

5.6.1 Hz Button

Pressing the Hz button toggles the Pulse Frequency Display mode. If Hz mode is on, the pulse frequency displays in the parameter edit region of the display. If Hz mode is off, the pulse frequency does not display. If the pulse frequency is >300 Hz, the displays shows a series of three dashes, followed by the letters, "Hz."

5.6.2 Zero Button

The Zero button is used to manually start a batch while in Statistics mode.


5.6.3 Power Switch and Backlight Toggle Button




The combination Power Switch and Backlight Toggle button serves a dual purpose. It sets power on/off to the meter, and toggles the backlight.

- When the meter is off, the power-on state is activated by pressing the button for one second.
- A press of the button for one second while in the power-on state sets the meter to OFF.
- When the meter is in the power-on state, the backlight state is toggled by pressing the button for less than one second.
- The backlight is always off when power is first applied to the meter.

5.6.4 Up and Down Arrows

 These buttons serve a dual purpose. When Edit mode is active, the buttons are used to adjust the currently-selected edit field. When Edit mode is not active, the buttons are used to select the measurement range.

5.6.5 Left and Right Arrows

 This button pair has a dual purpose. When the Edit mode is active, the buttons allow users to select the edit field of the currently-selected edit parameter. When the Edit mode is not active and Statistics mode is active, the buttons are used to select the statistical parameter of interest (MAX, MIN, MEAN). These buttons are nonfunctional when Edit mode and Statistics mode are not active.

5.7 Statistics Mode

FieldMaxII-P can be configured to display statistical data instead of instantaneous measurements. Statistical data is generated on a pulse-by-pulse basis.

For more information on selecting parameters while in Statistics mode, refer to 'Left and Right Arrows,' above.

The tuning needles and zone indicator bars are not present in Statistics mode.

5.7.1 Invalid Data

A statistical batch requires valid data for every data point in the batch. If a batch collection of data is in process and invalid data is measured, the batch is considered contaminated and the batch immediately ends with no batch result computed. The error that caused the contaminated batch is displayed

5.8 Measurement Display and Range Selection

The display update rate for numeric measurement is three times per second.

Measurement range is selected in decade steps. Range selection—shown in Table 5-2 (p. 30)—is dependent on the sensor characteristics, as well as user-determined measurement settings.

FieldMaxII-P uses the “3’s” Rule—a display formatting rule in which the display value is not allowed to exceed 3, 30, or 300, depending on where the decimal point falls, with the decimal point located in a fixed position, as determined by the current range. Typically if a reading exceeds the “3’s” limit, the instrument is over-ranged.

Over-ranging refers to a meter setup condition in which the sensor output signal is greater than the maximum allowable level for the selected range. An “OL” (overload) appearing on the display signifies an over-range condition. Over-ranging generates invalid data. Table 5-2 indicates the display format for different full-scale range settings.

Table 5-2. Full Scale Range Settings

Full Scale Measurement	Display Format
3 nJ	X.YYY nJ
30 nJ	XX.YY nJ
300 nJ	XXX.Y nJ
3 μJ	X.YYY μJ
30 μJ	XX.YY μJ
300 μJ	XXX.Y μJ
3 mJ	X.YYY mJ
30 mJ	XX.YY mJ
300 mJ	XXX.Y mJ
3 J	X.YYY J
30 J	XX.YY J
300 J	XXX.Y J

NOTICE

When area correction is enabled in Joules mode, a ‘/cm²’ will be appended to the units.s

The Range Hint annunciator displays the full-scale range value with the engineering prefix and units omitted. For example, when the range is 30 J, the Range Hint annunciator displays '30'. Note that it also displays '30' when the selected range is 30 mJ, 30 μ J, or 30 nJ. For more information about the Range Hint annunciator, refer to 'Range Hint' (p. 22).

5.8.1 Internal Triggering Mode

For greatest accuracy and repeatability, FieldMaxII-P must trigger reliably for each laser pulse. *Internal triggering* refers to extracting an artificial trigger from the incoming signal.

To successfully extract an internally-generated trigger, set the meter range so that the incoming signal is at least 5% of the full scale. Set the trigger level at least 2% above the noise level and at least 2% below the peak height.

For example, with a peak height of 300 mJ, set the meter to a range of 3 J. A peak will occur at 10% of full scale, well above the 5% level. The trigger level should not be set higher than 8%. If the noise level is 5% of full scale (approximately 150 mJ), the trigger level should not be set lower than 7%.

In the example shown in Figure 5-6, the internal trigger threshold has been set to 8% (shown as a dashed line). Pulse A will definitely not generate a reliable trigger. Pulse B may generate a trigger, but not reliably. Pulses C and D will definitely generate reliable triggers.

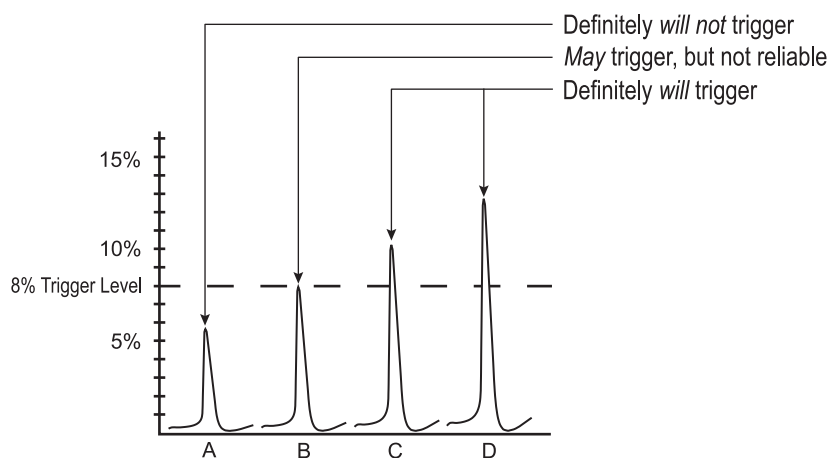


Figure 5-6. Internal Trigger Threshold

A full-scale signal on one range may not trigger on the next higher range unless the trigger level is set to less than 10%. For example, a near full-scale signal of 280 mJ on the 300 mJ range is less than 10% of full scale

on the next higher range (3 J), and would therefore not trigger the reading. To obtain reliable triggering in this instance, adjust the trigger level to less than 8%.

The trigger is synchronous with the leading edge of the pulse, but the actual peak is determined algorithmically by sampling the input signal near the trigger. From the trigger point forward, the algorithm searches for peaks and from the trigger point back, it searches for a baseline.

5.8.2 Hz Display Mode

If desired, pulse frequencies can be displayed when a pyroelectric sensor is attached to the meter. For more information, refer to 'Hz Button' (p. 28).

5.8.3 Zero the Sensor

The Zero button is used to manually re-zero the statistical batch.

6 Special Topics

This section gives more information on trigger states and modes of operation.

6.1 Trigger States and the Trigger Annunciator

Two segments are dedicated to the indication of the trigger state of the meter. These are the TRIG and the '?' segment that follows it.

There are two possible trigger states: 'triggered' and 'wait for trigger.' The triggered state indicates that the meter is capturing laser pulses. The wait for trigger state indicates an absence of pulse triggers. When a pulse is captured, the trigger state becomes active and the wait for trigger state becomes inactive. If no pulses are captured within 1.67 seconds since the last pulse was captured, the wait for trigger state becomes active and the triggered state becomes inactive. The TRIG annunciator reads 'TRIG' in the triggered state and 'TRIG?' in the wait for trigger state.

7 Host Interface

7.1 Overview

The host interface is intended for use with a National Instruments LabVIEW virtual instrument (VI) that accesses an ActiveX DLL called FieldMax2Lib.dll. The DLL handles communication and data transmission between the FieldMaxII and host applications written in LabVIEW.

A Getting Started introduction and LabVIEW examples are available at <https://www.coherent.com/resources>

8 Calibration and Warranty

8.1 Calibration

Coherent laser power and energy meters are precision instruments, capable of delivering very accurate measurements, as well as providing many years of useful service. To maintain this high level of performance, it is important to have measurement systems serviced and re-calibrated.

A large percentage of Coherent calibrations are performed within five business days, and expedited service is available, to minimize customer downtime.

8.1.1 Scope of Calibration

Calibration of a Coherent power and energy laser measurement product includes:

- Calibration to original uncertainty levels
- Minor repairs (see more information, below)
- Extended 12 month warranty, on eligible products
- Fast calibration turnaround time.

Minor repairs includes fixing manufacturer's defects, hardware updates, firmware, software updates, damaged connectors, and other small repairs. Detector element replacement due to laser damage and damage caused by negligent use is not covered – for customer caused damage, an additional repair service charge is applied.

This level of service results in an overall lower cost of ownership, with many owners realizing a lifetime warranty for their products.

8.1.2 Re-certify Once Per Year

To maintain this high level of performance, and to ensure compliance with your quality and ISO certification, it is important to have measurement systems serviced and re-certified once per year.

Extended use of laser power and energy meters and sensors, as well as environmental factors, can have an adverse effect on accuracy and also result in wear and/or damage to parts critical to optimum performance.

8.1.3 Coherent Calibration Facilities and Capabilities: ISO 17025 Accredited

Coherent calibration facilities contain the widest possible range of light sources from 193 nm to 10,600 nm, with powers ranging nanowatts to kilowatts.

Coherent Wilsonville, Oregon, and its satellite sites, are fully accredited to ISO/IEC 17025:2017 by ANAB - The ANSI National Accreditation Board. ANAB is a signatory of the International Accreditation Forum (IAF) and International Laboratory Accreditation Cooperation (ILAC) multilateral recognition arrangements. ANAB has signed the MRAs of the InterAmerican Accreditation Cooperation (IAAC) and the Asia Pacific Accreditation Cooperation (APAC). Accreditation to ISO/IEC 17025 is the formal recognition that a calibration laboratory is technically competent to carry out specific calibrations.

A detailed discussion of the Scope of Accreditation and the Technical Requirements of ISO 17025 Accreditation can be found on the Coherent ISO 17025:2017 Accreditation web page.

In addition, Coherent team delivers the industry's best service, with a knowledgeable and responsive staff, and rapid turnaround.

8.2 Limited Warranty

Coherent, Corp. (the "Company") warrants its laser power and energy meters and sensors products ("Products") to the original purchaser (the "Customer") that the product is free from defects in materials and workmanship and complies with all specifications, active at the time of purchase, for a period of twelve (12) months.

Coherent, Corp. will, at its option, repair or replace any product or component found to be defective during the warranty period. This warranty applies only to the original purchaser and is not transferable.

8.3 Warranty Limitations

The foregoing warranties shall not apply, and Coherent reserves the right to refuse warranty service, should malfunction or failure result from:

- Damage caused by improper installation, handling or use.
- Laser damage (including sensor elements damaged beyond repair).
- Failure to follow recommended maintenance procedures.

- Unauthorized product modification or repair.
- Operation outside the environmental specifications of the product.

Coherent assumes no liability for Customer-supplied material returned with Products for warranty service or recalibration.

THIS WARRANTY IS EXCLUSIVE IN LIEU OF ALL OTHER WARRANTIES WHETHER WRITTEN, ORAL, OR IMPLIED. COHERENT SPECIFICALLY DISCLAIMS THE IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE. IN NO EVENT SHALL THE COMPANY BE LIABLE FOR ANY INDIRECT, INCIDENTAL OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH ITS PRODUCTS.

9 Extended Warranty Program

Coherent, Corp. (the “Company”) offers original purchasers (the “Customer”) purchasing laser power and energy meters and sensors products (“Products”) an extended twelve (12) month Warranty program, which includes all parts and labor.

To qualify for this Warranty, a Customer must return the Product to the Company for recalibration and recertification.

- The Company will re-certify the Product, provide software upgrades, and perform any needed repairs, and recalibrate the Product, for a fixed service fee (as established by the Company from time to time and in effect at the time of service).
- If the product cannot be re-certified due to damage beyond repair, parts obsolescence, or other reasons, the Customer may be informed that an Extended Warranty program is not available for the Product.

If the Product fails and is returned to the Company within one year following the date of recalibration and recertification service, the Company will, at its option, repair or replace the Product or any component found to be defective. If the Product must be replaced and the Product is no longer available for sale, Coherent reserves the right to replace with an equivalent or better Product. This Warranty applies only to the original purchaser and is not transferable.

For information about calibration and recertification services, refer to ‘Calibration’ (p. 37).


Appendix A: Maintenance

A.1 Battery Replacement

FieldMaxII-P uses a rechargeable battery pack (standard), six 1.5V alkaline batteries, or a 90-to-260 VAC, 50/60 Hz AC adapter—refer to “AC Adapter” (p. 12) for more information. Figure A-1 (p. 44) illustrates how to replace the optional alkaline batteries.

Table A-1. Batteries Contained in this Product

Description	Type
7.2V rechargeable battery pack	NiMH
(optional) 1.5V AA	Alkaline

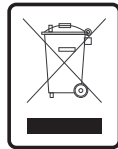

Dispose of batteries according to local regulations. Do not dispose as normal waste. Consult the local waste authorities for guidance.

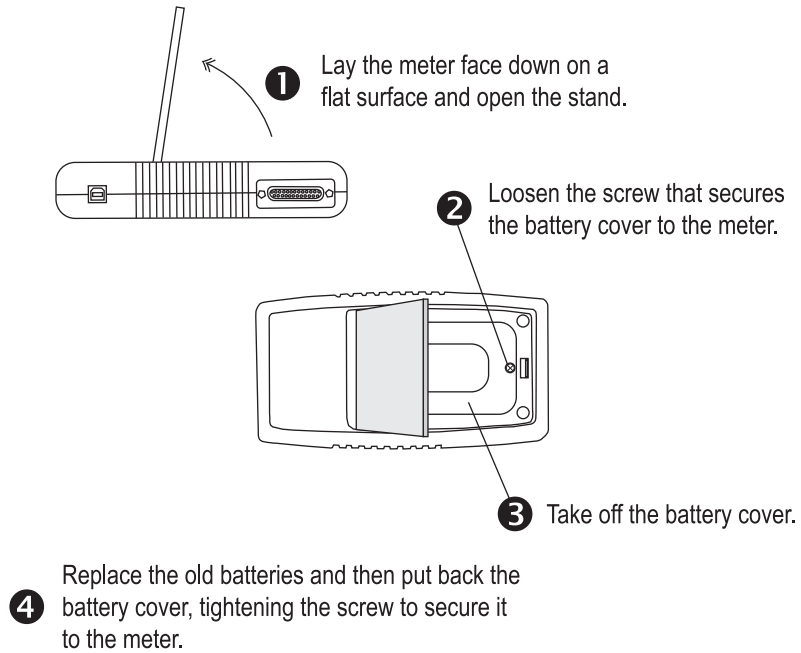


Figure A-1. Optional Alkaline Battery Replacement

Appendix A: Service & Support

This section provides information about technical support, service and shipping instructions.

A.1 Technical Support

Coherent provides telephone and web-based technical assistance as a service to its customers and assumes no liability for any injury or damage that can occur at the same time with such services.

Operation of any Coherent laser with any of its interlocks (or safety features) defeated is always at the operator's own risk. Under no circumstances do these support services affect the terms of any Warranty agreement between Coherent and the buyer.

Be prepared to provide the following information to the Product Support Engineer responding to the request:

- The unit model or part number
- Serial number
- A description of the problem
- Any corrective steps that have been attempted

A.1.1 Support in the USA and North America

Should any difficulties with the unit be experienced or for product or technical information, contact Coherent as follows:

- By email: LSMsupport@coherent.com
- Website: www.Coherent.com

For further assistance, please contact Coherent Technical Support:

- By phone: 1-(800)-343-4912 or 1-(503)-454-5700 outside the U.S.

Telephone coverage is available Monday through Friday (except U.S. holidays and company shutdowns). Inquiries received outside of normal office hours will be captured by our automatic answering system and calls will be quickly returned the next business day.

A.1.2 International Support

If located outside the U.S., visit www.Coherent.com for technical assistance, or contact a local Service Representative directly:

- Germany: +49-6071-968-0
- Japan: +813-5635-8680

On the Coherent website, contact information (telephone numbers and addresses) for Service Representatives worldwide is provided.

A.2 Obtain Service

To obtain product service under this warranty, Customer must notify the Company of the defect before the expiration of the warranty period and make suitable arrangements for the performance of service.

The Company shall, in its sole discretion, determine whether to perform warranty service at the Customer's facility, at the Company's facility, or at an authorized repair station.

If Customer is directed by the Company to ship the product to the Company or a repair station, Customer shall:

- Package the product (to protect from damage during shipping) as instructed in "Product Shipping Instructions" next.
- Ship it to the address specified by the Company, with shipping prepaid. back to Coherent in conjunction with recalibration and recertification.
- Coherent shall pay the cost of shipping the Product back to the Customer in conjunction with product failures within the first twelve (12) months of time of sale or during an extended 12-month warranty period.

A.3 Product Shipment Instructions

A Returned Material Authorization number (RMA) assigned by the Company must be provided on the outside of all shipping packages and containers. Items returned without an RMA number are subject to return to the sender. Detailed instructions to prepare a product for shipping are provided in the next section.

To prepare a product for shipping to Coherent:

1. Contact Customer Service for a Return Material Authorization number.
2. Attach a tag to the product that includes the name and address of the owner, the person to contact, the serial number, and the RMA number that was received from Coherent Customer Service. Pack this tag inside the box.
3. Wrap the product with polyethylene sheeting or equivalent material.
4. Using the original shipping and packaging materials, pack the product.
5. Seal the shipping carton with shipping tape or an industrial stapler.
6. Add the RMA number that was received from Coherent Customer Service to the shipping label on the outside of the box. Ship the product to one of the following addresses:

USA

Coherent Laser Measurement and Control Service Center
Attn: (your RMA number)
27650 SW 95th Avenue
Wilsonville, OR 97070

Europe

Coherent Shared Services B.V.
Dieselstr. 5 b
D-64807 Dieburg
Germany

Asia

Coherent, Inc. Japan
Atsugi Technical Center
Toda 1042-4
Atsugi, Kanagawa Prefecture
243-0023, Japan

Appendix B: Specifications

Table B-1, below, shows the specifications for the FieldMaxII-P.

Table B-1. Specifications

Parameter	Description
Electrical/Mechanical	
Analog Output	1, 2, or 5V full-scale (user-adjustable) 100 ohm source impedance
Battery Operating Time* (approx) * Without a backlight	<i>Rechargeable</i> <i>Alkaline</i> 8 hr. 24 hr.
Calibration Accuracy	± 1%
Digital Output	USB 1.1
Digital Tuning Needle	100 mS (tau) 20 Hz (update rate)
Power Requirements	<i>AC operation:</i> 90-to-260 VAC, 50/60 Hz <i>DC power input:</i> 12 VDC, 1.25A, center-positive <i>Battery operation:</i> rechargeable battery pack (standard) or six optional 1.5V AA alkaline batteries (battery specifications are listed earlier in this table under "Battery Operating Time")
Pyroelectric Input (maximum voltage input)	18V
Internal Trigger	2 to 20% full scale (selectable)
Linearity	± 1%
Measurement Resolution	± 0.1% of full scale
Environmental	
Altitude	4,500 m (operating) < 12,000 m (storage)

Table B-1. Specifications (Continued)

Parameter	Description
Relative Humidity	< 90% (5 to 40° C) (operating) < 95% (0 to 70° C) (storage)
Temperature	5 to 40° C (operating) -20 to 70° C (storage)
Ranges	
Area Parameter (entered as a diameter)	0.01 to 999.99 mm
Attenuation (Attenuation parameter)	0.01 to 999.99
Batch Size (Statistics parameter)	2 to 99,999 pulses
Measurement Range (full scale, sensor-dependent)	3 nJ to 300 kJ
Rep Rate	± 1 Hz (accuracy) 300 Hz (maximum) 1 Hz (resolution)
Window Size (Avg parameter)	2 to 99,999 pulses
Physical Characteristics	
Dimensions (h x w x d) (approx)	8 in. (20 cm) 4 in. (10 cm) 1.5 in. (4 cm)
Display	58 x 73 mm fixed-segment LCD with backlight
Weight (approx, including batteries)	1.1 lb. (0.5 kg)
Miscellaneous	
Regulations Met	CE

Index

A

- AC adapter 3-6
- Annunciator display functions 4-3
- Area button function 4-7
- ATTEN annunciator 4-4
- Atten button function 4-8
- AVG annunciator 4-4
- Avg button function 4-8

B

- Bad zero 4-9
- BATTERY annunciator 4-4
- Battery replacement 3-5
- Button functions
 - Area 4-7
 - Atten 4-8
 - Avg 4-8
 - Hz 4-9
 - Left and right arrows 4-9
 - Power switch and backlight toggle button 4-9
 - Setup/Local 4-6
 - Stat 4-6
 - Trig 4-8
 - Up and down arrows 4-9
 - Wave (λ) 4-7
 - Zero 4-9

C

- Calibration 7-1
- Coherent calibration facilities and capabilities 7-1
- Connector
 - Power jack 3-5
 - Sensor 3-5
 - USB 3-5

D

- Digital tuning feature in statistics mode 5-1
- Display functions
 - Annunciators 4-3
 - Fault displays 4-4
 - Invalid and not available data conditions 4-5
 - Measurement display and range selection 4-10

E

- Extended warranty 7-2

F

- Fault displays 4-4
- Front panel overview 3-2

H

- Hz
 - Button function 4-9
 - Display mode 4-12

I

- Internal triggering mode 4-11
- Invalid and not available data conditions 4-5
- Invalid data 4-10

L

- LCD display 3-3
- Left and right arrows 4-9
- Left side panel overview 3-5
- Limited warranty 7-2

M

- Measurement display and range selection display 4-10
- Meter modes
 - Hz display 4-12
 - Internal triggering 4-11

N

- Not available data 4-5

O

- Overview
 - AC adapter 3-6
 - Battery replacement 3-5
 - Front panel 3-2
 - Left side panel 3-5
 - Right side panel 3-4

P

- Power
 - Jack connector 3-5
 - Switch and backlight toggle button 4-9
- Product shipping instructions 7-4
- Pyroelectric-specific information
 - Hz display mode 4-12
 - Internal triggering mode 4-11

R

- Range hint annunciator 4-4
- Right side panel overview 3-4

S

- Sensor connector 3-5
- Service 7-3
- Setup parameters
 - Area 4-7
 - Atten 4-8
 - Avg 4-8
 - Setup/Local 4-6
 - Stat 4-6
 - TRIG 4-8
 - Wave 4-7
- Setup/Local button function 4-6
- Stat button function 4-6
- Statistics mode 4-10

FieldMaxII-P™

and digital tuning feature 5-1
and invalid data 4-10

T

Trig button function 4-8
Trigger states and the trigger annunciator 5-1
Tuning
 Mode 4-1
 Needles 4-2

U

Up and down arrows 4-9
USB connector 3-5




W

Warranty
 Extended 7-2
 Limitations 7-3
 Limited 7-2
Wave (lambda) button function 4-7

Z

Zero
 Bad 4-9
 Button function 4-9
Zone indicator bars 4-2

COHERENT

部件名称 Part Name	产品中有害物质的名称及含量						  
	有害物质 Hazardous Substances						
	铅 (Pb)	汞 (Hg)	镉 (Cd)	六价铬 (Cr(VI))	多溴联苯 (PBB)	多溴二苯醚 (PBDE)	
印刷电路板组装 Printed Circuit Board Assembly	X	O	O	O	O	O	
电源 Power Supply	X	O	O	O	O	O	
电源线 Power Cord	X	O	O	O	O	O	

本表格依据 SJ/T 11364 的规定编制
 O: 表示该有害物质在该部件所有均质材料中的含量均在 GB/T 26572 规定的限量要求以下。
 X: 表示该有害物质至少在该部件的某一均质材料中的含量超出 GB/T 26572 规定的限量要求。

Download software and manuals at <https://www.Coherent.com/>



FieldMaxII-P™ User Manual
 ©Coherent Corp., 2024 (RoHS), printed in the USA
 Part No. 1086242 Rev. AH